


Esco Sample Protection and Remote Monitoring Solutions *Sample monitoring made easy*

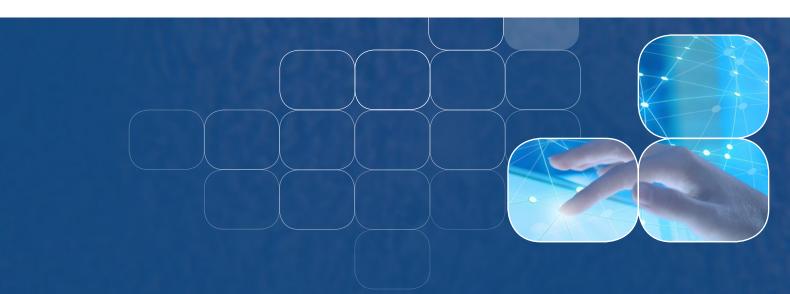

- **Voyager**® (Remote Monitoring, Programming, Datalogging Software)
- **iRMAP**TM (Intelligent Remote Monitoring Application Protocol)
- **Voyager® Box** (Web Monitoring and Datalogging System)
- **PROtect** (A Completely Independent and Redundant Sample Monitoring System)
- SampleTracker (Biosample Inventory Management System)



Esco Sample Protection and Remote Monitoring Solutions

Table of Contents

Corporate Profile	3
Products and Applications	4
Introduction	5
Voyager® Remote Monitoring, Programming, Datalogging Software	6
Voyager® Key Features	7
Voyager® Interface with Esco Equipment	8
Voyager® Installation and System Requirements	9
iRMAP™ Intelligent Remote Monitoring Applications Protocol	. 10
iRMAP™ Key Features	11
Voyager® Box	13
Voyager® Box Features	14
Voyager® Box Installation and System Requirements	15
PROtect A Complete Independent and Redundant Sample Monitoring System	16
PROtect Key Features	. 17
PROtect System Components and Requirements	. 19
PROtect General Specification	20
PROtect Options and Accessories	21
SampleTracker	24
SampleTracker Key Features	25
Esco After sales Service	27
Esco Global Network	28





Welcome to Esco

Esco's Vision is to provide enabling technologies for scientific discoveries to make human lives healthier and safer.









The Esco Group of Companies is committed to deliver innovative solutions for the clinical, life sciences, research, industrial, laboratory, pharmaceutical, and IVF community. With the most extensive product line in the industry, Esco have passed a number of international standards and certifications, and is operating under ISO 90001, ISO 14001, and ISO 13485. Esco represents innovation and forward-thinking designs, that are of the highest standard quality since 1978.

Availability and Accessibility. Esco has headquarters in Singapore, Indonesia, and Philippines, with manufacturing facilities are located in Asia and Europe. Research and Development (R&D) is conducted worldwide spanning the US, Europe and Asia. Sales, services and marketing subsidiaries are located in 42 major markets including US, UK, Japan, China and India. Esco regional distribution centers are located in Singapore, Malaysia, Thailand, Vietnam, Myanmar, Indonesia, Philippines, Bangladesh, Hong Kong, Taiwan, South Korea, China, Japan, India, UAE, Central and South Africa, Denmark, Germany, Italy, Lithuania, Russia, United Kingdom, and USA. Because of our worldwide presence, you can be sure that Esco is within your reach.

High Quality, Reliable, and Dependable. Esco products are of high quality, reliable, and dependable; assuring customers of research accuracy. Cross functional teams from Esco Production, R&D, Quality Assurance, and Senior Management, are regularly assembled to review and implement areas for improvement.

Esco Cares for Your Safety. Esco focuses on providing safety not just for your samples but also for you and the environment.

Esco Cares for Your Comfort. Building ergonomic designs and reducing noise levels of the units ensures comfort for our users.

Esco Cares for the Environment. One in every four of Esco's employees is involved in R&D and a number of them evaluate new components and/or designs to produce energy efficient equipment. Being GREEN is more than just modifying parts used to produce a new energy efficient technology, it is also embodied in the every aspect of the company.

Customer Service and Support. Our service does not stop once purchase has been done. Esco gives on-time customer service and offers end-user seminars, service training, preventive maintenance, and provides educational materials and informative videos.

As Esco takes the opportunity to respond to the world's needs, we aim not only to contribute in the advancement of scientific discoveries but also in making the world a safer, healthier, and better place to live in.

Products and Application

Life Sciences Laboratory Equipment

Sample Preparation

- Class I Biological Safety Cabinets
- Class II Microbiological Safety Cabinets
- Class II Type A2 Biological Safety Cabinets
- Class II Type B1 Biological Safety Cabinets
- Class II Type B2 Biological Safety Cabinets
- Class III Biological Safety Cabinets
- Horizontal Laminar Flow Clean Benches
- Vertical Laminar Flow Clean Benches
- Laboratory Animal Research Workstations
- Laboratory Centrifuges

Sample Cultivation

- CO, Incubators, Direct Heat Air-Jacketed
- CO₂ Incubators with Cooling System
- CO₂ Incubators with Stainless Steel Exterior
- Laboratory Shakers

Sample Handling and Analysis

PCR Thermal Cyclers

• Conventional Thermal Cyclers

PCR Sample Handling

- Microplate Shakers
- PCR Cabinets

Sample Storage & Sample Protection Solutions

- Ultra-low Temperature Freezers
- Lab Refrigerators and Freezers
- Sample Database Management Software
- Intelligent Remote Monitoring Application Protocol
- Remote Monitoring, Datalogging, Programming Software
- Wireless Monitoring System

Chemical Research

- Ductless Fume Hoods
- Laboratory Fume Hoods
- Fume Hood Airflow Monitors
- Exhaust Blowers
- Powder Weighing Balance Enclosures
- Filtered Storage Cabinet

General Equipment

Laboratory Thermostatic Products

- Laboratory Oven
- · Laboratory Incubator
- · Refrigerated Incubator
- Natural Convection Incubator

Forensic Sciences

• Evidence Drying Cabinet

Medical / IVF Equipment

Controlled Embryo Handling

- Fertilisafe™ ART Workstation
- AVT-I Anti Vibration Table
- Semi Closed Environment IVF

Safe Embryo Culture

- MIRI® Multi room Benchtop Incubator
- CelCulture® CO2 Incubator
- Mini MIRI[®] Humidified Benchtop Incubator

Innovative Time Lapse Imaging

• MIRI® Time-lapse Incubator

Accurate Quality Control

- MIRI® GA Gas and Temperature Validation Unit
- MIRI® GA Mini Gas Validation Unit

Unique Consumables

CultureCoin

Esco Pharma Products

Airflow Containment Products

- Pharmacon® Downflow Booths
- Ceiling Laminar Airflow Units
- Laminar Flow Horizontal/Vertical Trolley
- Enterprise™ Laminar Flow Straddle Unites
- Cytoculture[™] Cytotoxic Safety Cabinets

Isolation Containment

- Aseptic Containment Isolator (ACTI)
- Weighing and Dispensing Containment Isolator (WDCI)
- General Processing Platform Isolator (GPPI)
- Containment Barrier Isolator (CBI)
- Turbulent Flow Aseptic (Grade A) Isolator (TFAI)
- Isoclean® Healthcare Platform Isolator (HPI)
- Streamline® Compounding Isolators (SCI)
- Technetium Dispensing Isolators
- Blood Cell Labeling Isolators
- Open and Closed Restricted Barrier Access Systems (RABS)

Cross Contamination Facility Integrated Barrier

- BioPass™ Pass Through
- Infinity® Air Shower Pass Box
- Cleanroom Air Shower
- Infinity® Cleanroom Transfer Hatch
- Infinity[®] Pass Box
- Soft capsule® Soft Wall Cleanroom
- Dynamic Passboxes and Dynamic Floor Laminar
- Laminar Flow Storage Cabinet

Ventilation Containment

• Ventilated Balance Enclosure

VacciXcell Products

Bioreactors and Fermenters

- CelCradle™
- TideXcellTM
- VacciXcell™ Hybrid Bioreactor

Cell Culture Monitoring, Media and Consumables

- Super Plus™
- Plus™ Vero
- Plus™ MDCK
- Plus™ MDCK II
- BioNOC™ II macrocarriers
- GlucCell™ Glucose Monitoring System
- CVD Kit

Filling Line Equipment

- Filling Line Isolators
- cRabs (close restricted access barriers)
- oRabs (open restricted access barriers)

Integrated Solutions

- Cell Processing Isolator • Cell Processing Center

TaPestle Rx Products and Services PRODUCTS

Pharmacy Automation and Compounding Supply

- Compounding Pharmacy Isolators (SCI, HPI, CBI, GPPI)
- Safety Cabinets and Enclosures (Class II BSC, VBE, LFC)
- Radiopharmacy Hoods and Isolators
- Aseptic Filling Systems

Healthcare and Laboratory Construction Components

- Prefabricated Walls (Airecell®)
- Prefabricated Containerized Facility (Prefab™)
- Series Ceiling Systems
- Hygienic/Hermetic Door Systems
- Surgical Scrub Sinks
- Vinyl Tiles and Epoxy
- Laboratory Fit-outs
 - Worktops - Frames
 - Specialty Storage cabinets
 - Service Spines & Reagent Shelving


SERVICES

- Conceptualization
- Planning
- Procurement Installation
- **FACILITY DESIGNS**
- Process Architecture
- Biocontainment/Biosafety
- Pharmacy Compounding/Nuclear Medicine • Cleanroom, Vaccine and Cell Processing
- Laboratory
- Containerized Facility
- ART/IVF
- Cold Chain

Every precaution has to be taken to protect a sample's integrity.

Be assured with the guaranteed quality of performance and service Esco can provide.

Esco's Sample Protection and Remote Monitoring Solutions provides an extensive solution to prevent loss and spoilage of unique samples and products.

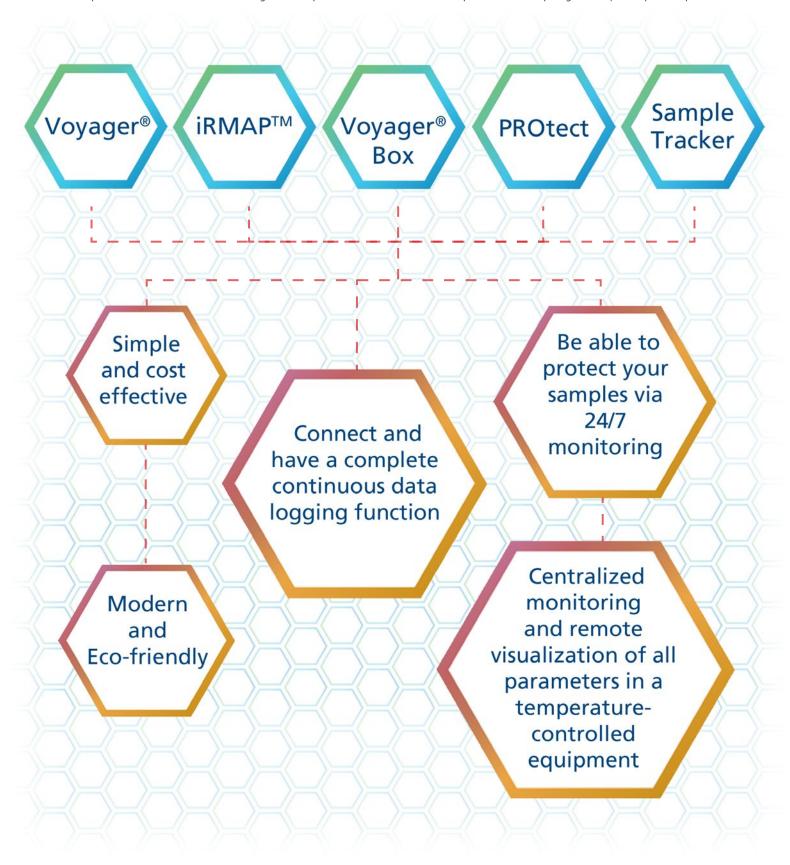



Voyager®

Remote Monitoring, Programming, Datalogging Software

Esco Voyager® is a PC-based software package developed for the remote monitoring, datalogging, and programming/ device configuration of Esco controlled environment laboratory equipment. The equipment has the capability to push updates to Esco equipment connected to software. In any event of equipment failure, the software is able to send out remote alarm notification through e-mail or SMS with a modem configured. It is also compliant to 21 CFR Part 11 standard for traceability.



$\textbf{Compatible Esco Equipment} \ (\textit{Clockwise from top}):$

Lexicon® Ultra-low Temperature Freezer (UUS) Isotherm® Refrigerated Incubator (IFC) Isotherm® Forced Convection Incubator (IFA) CelCulture® CO, Incubator (CCL)




KEY FEATURES




Remote Monitoring

- Automatic, continuous monitoring of device parameters (Temperature, CO₂, RH, etc.) to validate instrument performance and assay criteria
- Automatic e-mailing of log data at user-defined intervals

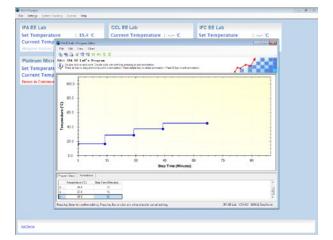

Remote Programming and Equipment Configuration

- Development of programs using a graphical interface, for subsequent download to device memory
- Voyager® software can download user programs to multiple Esco instruments easily insuring consistent setup on multiple devices.
- Configure devices remotely (measurement interval, temperature set point, low temperature alarm point and high temperature alarm point)
- This window also allows the user to choose the method of when and which data can be viewed (spreadsheet, email, web page).

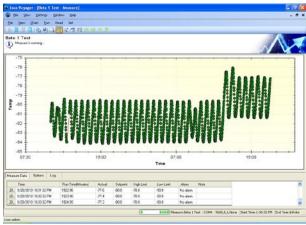

Datalogging and Graph Functions

- Easy viewing and graphing multiple device parameters remotely in real-time
- Saving and exporting log data in various formats from multiple instruments
- Download of data logs from equipment memory (only for devices with built-in data logging memory, i.e. CO₂ Incubators)

Alarm

- Alarm is activated if a user-defined parameter limit is exceeded.
- Automatic email alerts
- Documentation of all alarms with time and date stamp





Voyager® interfaces with individual Esco equipment over various available communication ports with the aid of EscoBUS communications protocol.


Compatible Equipment

- Lexicon II® Ultra-Low Temperature Freezer (Silver, Gold, and Platinum)
- CelCulture® CO₂ Incubator (CCL)
- CelMate® CO₂ Incubator (CLM)
- Isotherm® Forced Convection Laboratory Oven

- Isotherm® Natural Convection Laboratory Incubator (INA)
- Isotherm® Forced Convection Laboratory Incubator (IFA)
- Isotherm® Refrigerated Incubator (IFC)



Installation Requirements

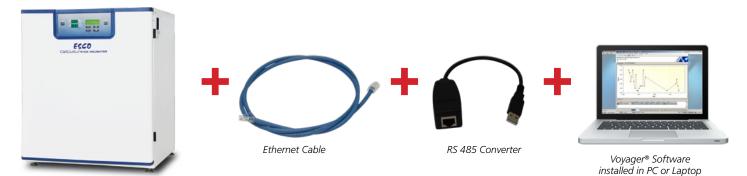

Compatible Equipment

- Industrial strength, industry-standard RS485
- EscoBUS communications protocol
- RJ45 ports at rear of each piece of equipment at the rear
- RJ45 to USB interface, plugs into PC USB port
- RJ45 to RS232 interface, plugs into PC DB9/Serial Port
- Each device on the network is assigned with a unique EscoBUS address
- Up to 4 devices of CO₂ Incubator may be daisy-chained to a single PC
- Up to 32 combined devices may be daisy-chained to a single PC
- The maximum distance between equipment is 100 m (328 ft), while the maximum distance between the PC to the last equipment is 1000 m (3280 ft).

Minimum System Requirements

Operation system: 32-bit or 64-bit Windows 2000, 2003 Server, XP, 7 Professional, and 10.

Memory: 4 GB Hard disk space: 10 GB Interface: USB/RS232/ RS485

Prerequisite Software: .NET Framework 2.0 (included)

Windows Installer 3.1 (included)

^{*}Internet connection required for email alert feature.

ORDERING INFORMATION		
Software	Item Code	Description
Voyager® and iRMAP™	5250001	Voyager® Software Kit with iRMAP™ subscription for 2 years







$iRMAP^{TM}$

(Intelligent Remote Monitoring Application Protocol)

24/7 Sample Monitoring

Intelligent Remote Monitoring Application Protocol (iRMAP TM) is a software system for remote monitoring of Voyager $^{\circ}$ -compatible Esco devices. A user purchases Voyager $^{\circ}$, installs it on a PC and connects it up to the network of Esco equipment. The monitoring data is encrypted and transmitted to an Esco server. User will be able to monitor the freezers/ cabinets 24 hours a day, 7 days a week from anywhere in the world.

The iRMAP, in conjunction with the Predictive Failure Monitoring System can also monitor the health of the Esco unit and proactively dispatch service to correct pending and future issues that could threaten customer samples. iRMAP™ can be accessed anytime, anywhere using a web browser/ iOS/ android mobile application.



KEY FEATURES

Sample monitoring anytime, anywhere

- 24/7 sample monitoring via internet browser and mobile iOS/ android application
- In the event that your samples are at risk, user will be consistently notified until situation is resolved.
- Every connected Esco unit's health will be monitored and service will be proactively dispatched to correct pending and future issues that could threaten customer samples.










Accessibility

• Users can access Voyager data logs anytime and at any location using either web browser applications, iOS, and android mobile devices.

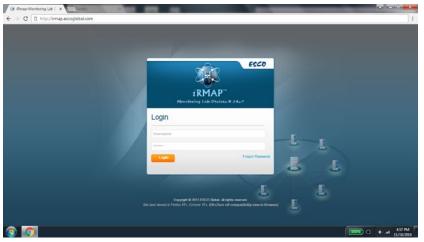










iOS Phones (Beta)



Alarm


- Alarm is activated if a user-defined parameter limit has exceeded.
- On-screen display of alarm notification and email prompts
- Documentation of all alarms with time and date stamp

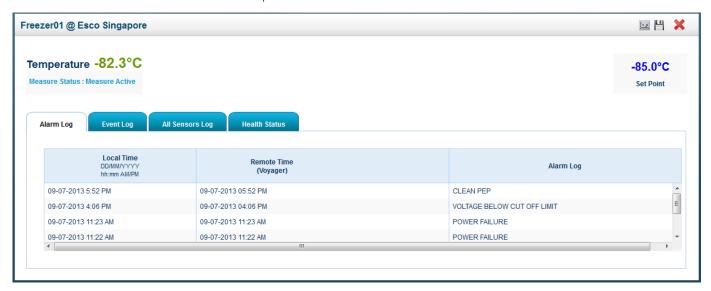

Data Log History Display and Graph Functions

- Remote viewing of data log history from multiple devices
- Actual data graph, event and alarm logs history
- Easy export of data as Excel csv file



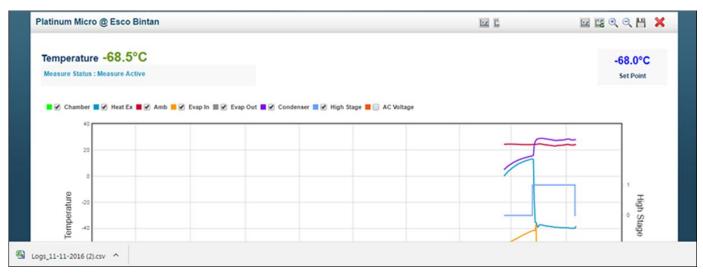

CLASS. WORLDWIDE.

Voyager® Box

System Overview

Esco Voyager Box is a web-based application developed for remote monitoring, and datalogging of Esco controlled environment laboratory equipment. The device will enables you to monitor your important equipments wirelessly, wireless Local Area Network (LAN) is needed to access the monitored unit. Each monitored parameters can be viewed using web browser of PC, laptop or smartphone. In any event of equipment failure, the application is able to send out remote alarm notification through SMS or e-mail.




Softwares and Applications

Compatible Equipment

- Lexicon II® Ultra-Low Temperature Freezer (Silver, Gold, and Platinum)
- CelCulture® CO2 Incubator (CCL)
- CelMate® CO2 Incubator (CLM)
- Isotherm® Forced Convection Laboratory Oven
- Isotherm® Natural Convection Laboratory Incubator (INA)
- Isotherm® Forced Convection Laboratory Incubator (IFA)
- Isotherm® Refrigerated Incubator (IFC)



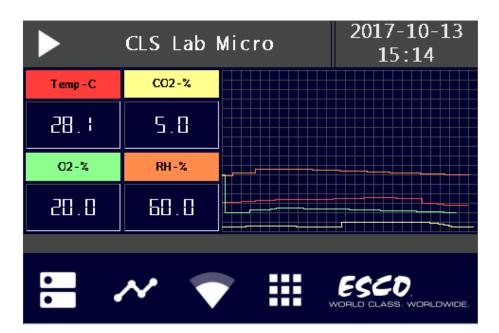

Home Menu

- Touch screen 3,5" display
- Display all monitored parameters
- Access all menu, Wifi connection, add device

Device		
CLS Lab Micro		
CCL Lab Electro		
IF A Lab Biology		
IFC Lab		
ULTF Lab		Ť
← ±	*	×


Add Devices

- List monitored device
- Configure measurement
- Add, edit, and delete monitored device

Measur ement			
Time	Log		
2017-10-13 08:05	Temp: 28.1/CO2:5.0/O2:20.0/RH:60		
2017-10-13 08:07	Temp: 27.8/CO2:5.0/O2:20.1/RH:61		
2017-10-13 08:09	Temp: 28.1/CO2:5.0/O2:20.1/RH:61		
2017-10-13 08:11	Temp: 28.0/CO2: 5.0/O2: 20.0/RH: 60	Ť	
←	• •		

Log Measurement

- Measurement history
- Up to 365 days measurement
- Save logging data with USB

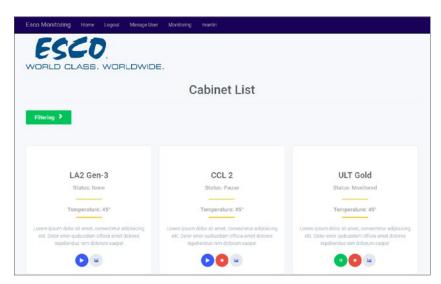


Web Interface

- Access all the monitored device in web browser
- Configure the measurement settings as simply as in VoyagerBox device



Installation Requirements

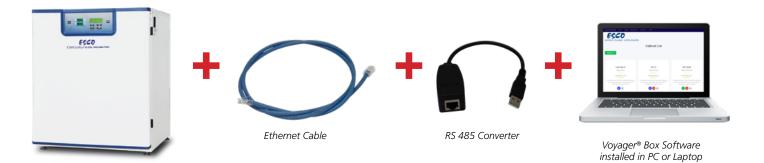

Minimum System Requirements

Operation system: 32-bit or 64-bit Windows 2000, XP, 7

Memory: 512 MB Hard disk space: 1 GB Interface: Wifi Connection

Compatible Equipment





PROtect

A Completely Independent and Redundant Sample Monitoring System


Esco PROtect monitoring system includes radio frequency recorders which can be connected to any brands of freezer, cold storage system and temperature controlled equipment. These recorders automatically send data to a central server to monitor temperature and other parameters and send real-time alert to users in case of severe failure. Esco PROtect complies with quality assurance procedures that is ISO 17025, ISO 15189 norm, GMP and audit trail traceability for conformance to 21 CFR Part 11.

Esco PROtect Wireless Monitoring System can be installed in the following industries/ sectors:

- Hospitals
- Pharmaceutical Industry
- Blood Bank Centers
- Food Safety Laboratories

- Animal Testing Laboratories
- Drugstores
- Research Centers
- Bio-Banking and Repositories

- Private Clinical Laboratories
- Food Industries
- Universities/Academia
- Cosmetic Industry

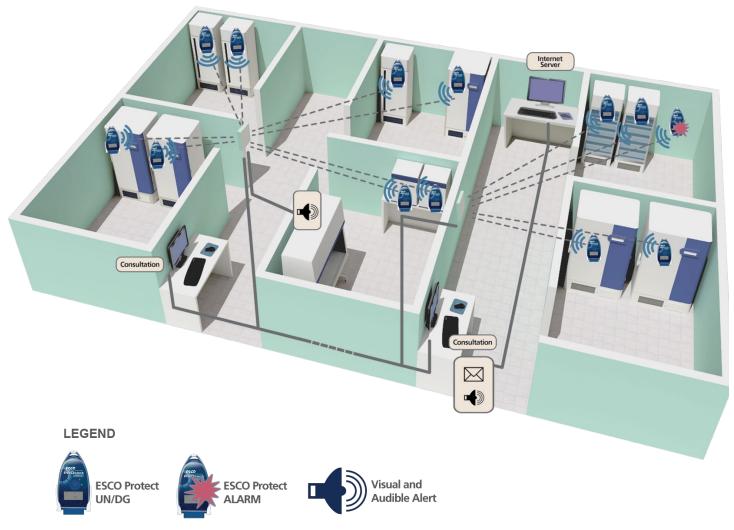



ESCO Protect EMD-1



Email Alert _

__ Ethernet network ______ Radio Communication (200-300m)

KEY FEATURES





Real Time Monitoring System

- Wireless Technology that protects your samples 24/7 with continuous data logging function and immediate alarm notifications
- Centralize monitoring for all your equipment (any brands and any models):
 - Temperature for refrigerators, freezers, cold chambers, cryogenic tanks and -80°C freezer
 - Temperature and CO₂ for ovens/incubators
 - Temperature and humidity for climatic chamber, ambient area
- Measurement and recording of physical parameters such as temperature, CO₂, and humidity
- Direct reading on the display
- Radio frequency has longer battery life and better transmission distance, as compared to Wi-Fi
- In using the radio frequency, repeaters can be installed to extend the distance which cannot be easily done by some Wi-Fi system

Simple and Cost Effectiveness

- Fast Installation, commissioning and maintenance
- Can centrally monitor integrity of samples even during a power outage or electronic failure of refrigerator, freezer, incubator, oven, etc
- Accurate and efficient compared to manual data readings
- Reduce labor cost for manual data readings. Manual data readings are less accurate, less reliable and time consuming. Saves you time and money.
- Readily accessible thru the display or monitoring software thru PC
- Multi-site installation
- Interchangeable digital probes with embedded memory to facilitate maintenance procedures.
- Modern and eco-friendly

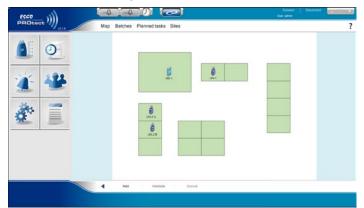

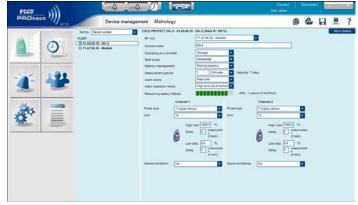

Traceability and Security - Quality Assurance

 Comply with quality assurance procedures: ISO 17025 and ISO 15189 norms, Good laboratory practices and local regulations recommend the use of tools to guarantee the reliability of your temperature controlled equipment (ultra-low temperature freezers, lab freezers, lab refrigerators, lab ovens, lab incubators, refrigerated incubators, CO₂ incubators, etc).





- Audit trail traceability for conformance to 21 CFR part 11
- Local regulations or recommendations within Quality Management Programs





Alarms

- Real time alarm management
- Warns if problem occurs via audible and visual alarm devices
- Remote alarm notification via SMS/ E-mail









Remote Monitoring and Datalogging

- Remote visualization of data
- Downloaded database is presented in easy-to-understand reports or graphs
- PROtect software enables to visualize recorders located on the installation map and to manage your alarm

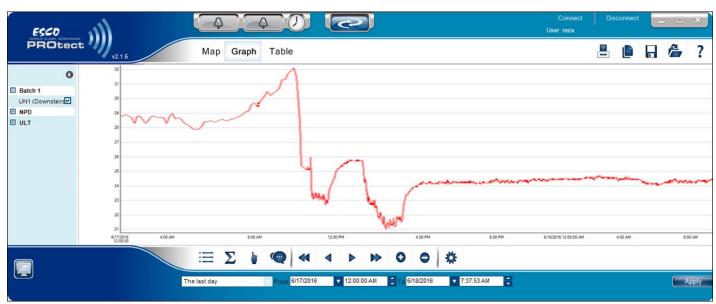

System Components



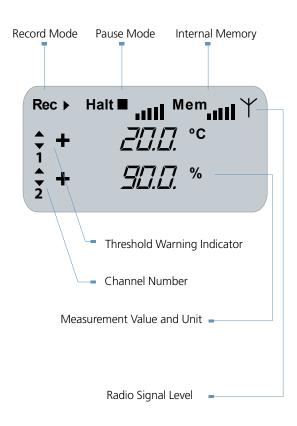

Potential Users of Esco PROtect Wireless Monitoring System

- Hospitals
- Pharmaceutical Industry
- Blood Bank Centers
- Food Safety Laboratories
- Animal Testing Laboratories
- Drugstores
- Hospitals
- Pharmaceutical Industry

- Private Clinical Laboratories
- Food Industries
- Universities/Academia
- Cosmetic Industry

Minimum System Requirements

PROtect Server:

Computer Architecture: 32 bit or 64 bit

OS: Windows 2000, XP, Vista, Windows 7 Professional and Windows 2003-2012 Server

RAM: 8 GB

HD: 250 GB capacity

Video Card: VGA compatible 1280 X 1024 32 bit

PROtect Client:

Computer Architecture: 32 bit or 64 bit

OS: Windows 2000, XP, Vista, Windows 7 Professional, Windows 2003-2012 Server and Windows 10

RAM: 4 GB

HD: 20 GB capacity

Video Card: VGA compatible 1280 X 1024 32 bit

PROtect can be used on the following regions:

- Europe 868 Mhz
- Africa 868 Mhz
- Asia 868 Mhz
- India 868 MhzANZ 902 Mhz
- North America 902 Mhz
- South America- 868 Mhz/902 Mhz
- Middle East 868 Mhz
- * Currently not applicable for South Korea and Japan region





GENERAL SPECIFICATION	Esco PROtect UN	Esco PROtect DG	
Measurement Range	From -200 °C to + 400 °C / 4-20mA / 0-1V /dry contact	T:-40 °C to +80 °C / tH:-30 °C to +70 °C from 0 to 100% RH non-condensing	
Channel Number	1 or 2	1 or 2	
Resolution	0,1℃	0,1℃	
Operating Conditions	From -30°C to +70°C	From -30°C to +70°C	
Power Supply	Lithium battery	Lithium battery	
Battery Lifetime	2 years - Replaceable	2 years - Replaceable	
Type of Input	External (Pt 100/Pt 1000/CO ₂ Probe/ 4-20mA/0-20mA/0-1V/dry contact)	External (temperature, relative humidity and dry contact)	
Recording Interval	1 second to 90 minutes	1 second to 90 minutes	
Internal Memory	20 000 measurements	20 000 measurements	
Radio Range	300 m	300 m	
Radio Bande	868 Mhz or 902 Mhz ISM	868 Mhz or 902 Mhz ISM	
Power	25 mW	25 mW	
Digital Display	Yes	Yes	
Dimension	123 x 69 x 35 mm	123 x 69 x 35 mm	
Protection Index	IP65	IP65	
Conformities	EN 12 830, ROHS, CE, FCC, ETS 300-220	EN 12 830, ROHS, CE, FCC, ETS 300-220	
Part number	868 Mhz UN-1: 1081050 UN-2: 1081052 902 Mhz UN-1: 1081051 UN-2: 1081053	868 Mhz DG-1: 1081054 DG-2: 1081056 902 Mhz DG-1: 1081055 DG-2: 1081057	
Included	User manual, wall-mounting bracket and cable input protection	User manual, wall-mounting bracket and cable input protection	



Ordering Information

Modems, Relays and Alarm System			
Product Image	Item Codes	Description	Features/Specification
PROTECT	1081050	Esco PROtect UN-1 868 Mhz band	• Single Channel RF Transmitter, -200°C to + 400°C; 4-20mA; 0.1V;
	1081051	Esco PROtect UN-1 902 Mhz band	Dry Contact
	1081052	Esco PROtect UN-2 868 Mhz band	• Dual Channel RF Transmitter, -200°C to + 400°C; 4-20mA; 0.1V ;
	1081053	Esco PROtect UN-2 902 Mhz band	Dry Contact
	1081054	Esco PROtect DG-1 868 Mhz band	• Single Channel RF Transmitter, -40°C to +80°C / tH:-30°C to +70°C
PROTECT COST	1081055	Esco PROtect DG-1 902 Mhz band	from 0 to 100% RH non-condensing
	1081056	Esco PROtect DG-2 868 Mhz band	• Dual Channel RF Transmitter, -40°C to +80°C / tH : -30°C to +70°C
	1081057	Esco PROtect DG-2 902 Mhz band	from 0 to 100% RH non-condensing
FICE	1081058	Esco PROtect MD-1 868 Mhz band	 Connects to your PC, it enables communication with Esco PROtect UN/DG Communicates with an unlimited number of recorders. Remote alarm driven by integrated dry contact output (vocal telephonic transmitter, buzzer, light).
	1081059	Esco PROtect MD-1 902 Mhz band	Software, PC or power supply failure alarm. Can be used with or without Esco PROtect Relay.
E5CO	1081060	Esco PROtect EMD-1 868 Mhz band	• Enables Ethernet connection to maximize distances between recorders
FSC0 PROtect	1081061	Esco PROtect EMD-1 902 Mhz band	and PC servers
FROTEST	1081062	Esco PROtect RL-1 868 Mhz band	 Enables direct RF communication to maximize distances between recorders and PC servers. Works with external power supply and supplied with a rechargeable battery.
	1081063	Esco PROtect RL-1 902 Mhz band	Includes two power ranges: a) 25 mW to communicate with recorders b) 500 mW to communicate with Modem
FROTOGE	1081064	Esco PROtect AM-1 868 Mhz band	 Provides real time warning if a problem occurs Integrated sound and visual alarm. Works with an external power supply and supplied with a backup battery.
	1081065	Esco PROtect AM-1 902 Mhz band	 Push button for local alarm acknowledgement. Relay output to trigger remote alarm (vocal telephonic transmitter) Warns Esco PROtect software when the battery is low.

Ordering Information

Options and Accessories			
Product Image	Item Codes	Description	Features/Specification
RIC SAME	1081066	PROtect Test-1 868 MHz	Use as a recorder to test connection between modem and recorder with
	1081067	PROtect Test-1 902 MHz	actual percentage signal
PICA PRODUCT	1081068	PROtect Modem Test-1 868 MHz	Use as a modem to test connection between modem and recorder with
(as No. 72)	1081069	PROtect Modem Test-1 902 MHz	actual percentage signal
	1081070	PROtect Software and Dongle	 Remote visualization of data. Downloaded database is presented in easy to understand reports or graphs. PROtect software enables to visualize recorders located on the installation map and to manage your alarm. Compliant with FDA (CFR 21 part 11) Note: PROtect software and dongle needs to be installed even if one (1) PROtect recorder (UN or DG) is connected to the equipment.
	1081071	Audible and Visual Alarm Device	 Warns if an alarm occurs. Works with an external power supply. Connects directly to PROtect Modem or to PROtect Alarm. Audible and visual Alarm.
	1081072	PT100 Probe Class B	
	1081288	PT100 Probe Class A	Measurement range: -200°C to + 260°CType of cable: 3 wires (3m)
	1081289	PT100 Probe Class 1/3 DIN	,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,,
	1081073	Digital Temperature Probe	 Type of probe: Dallas Measurement range: -40°C to + 80°C (-45°C max) accuracy: ± 0.3°C Type of cable: flat 3m cable Embedded memory for gauging parameters storage
	1081075	Battery for Recorders	• Lithium Battery for 10 pcs
	1081170	Digital Temperature / Humidity Probe	• Measurement range : -40 +80°C /0 100%HR
• • •	1081169	Digital Temperature/ Humidity Probe with 3m Extension Lead	 Type of probe for T°: Pt 100 Class A Class A Accuracy: ±2%HR from 0 to 90%HR ±3%HR from 90 to 100%HR
	1081293	Digital Temperature/ Humidity Probe with 12m Extension Lead	Power: delivered by the SPYRF recorder
	1081166	Digital (Low) Temperature	 Type of probe: Low Temp Stainless Steel probe Type of cable: 3m Teflon Cable Measurement range from 0°C to -200°C accuracy: ± 0.2°C Immersion of cable: continuous immersion in water and glycolic solutions and punctual immersion in alcoholic solutions"
	1081080	Universal Adaptor Cable for CO ₂ and DP sensors	• (5m) with connector IP67 for PROtect UN and DG.

Ordering Information

Options and Accessories			
Product Image	Item Codes	Description	Features/Specification
	1081165	Extension Lead for Probe	• 3m Extension Lead for Digital Temperature Probe
	1081262	SMS Modem	 Warns if any alarm occurs by sending SMS to receiver's number Needs a SIM card Supports multiple receiver's number USB Connection
	1081076	PROtect CO ₂ Probe	 Measurement range: 0 – 20% Output: 4-20 mA Mounting: Wall mounting bracket supplied Supply: 100-240 VAC
	5072029	PROtect Differential Pressure Probe	 Measurement Range: -100 Pa to +100 Pa Output: 4 - 20 mA Supply: 100-240 VAC (include Power Supply 1081263)
	1081078	Power Supply for Modem	• Power Supply for MD-1, EMD-1, AM-1
7	1081079	Power Supply for Relay	• Power suppy for RL-1
	1081167	PROtect Door Contact (With Sliding Doors)	• Use to sense door contact that door is open or close
	1081168	PROtect Door Contact (Non-Sliding Doors)	Use to sense door contact that door is open or close
COM Auto-males	5072009	PROtect GSM Auto Dialer set	 Warns if any alarm occurs by call receiver number (auto dialer) Needs a SIM card Supports until three (3) receiver numbers Supply: 100-240 VAC (include Power Supply 1081291)






SampleTracker


System Overview

SampleTracker is a comprehensive Bio sample Inventory Management solution. The Software tracks the sample management activities throughout its life cycle. The software scope covers all the features and functionalities required for the Sample Inventory Management and Tracking for all sizes of laboratories, managing different types of samples and having various types of storages. Flexibility in configuration of laboratories and sample attributes is one of the unique characteristics of SampleTracker.

SampleTracker automatically generates a unique identification number for each sample, which makes it easy for users to trace the sample as and when required

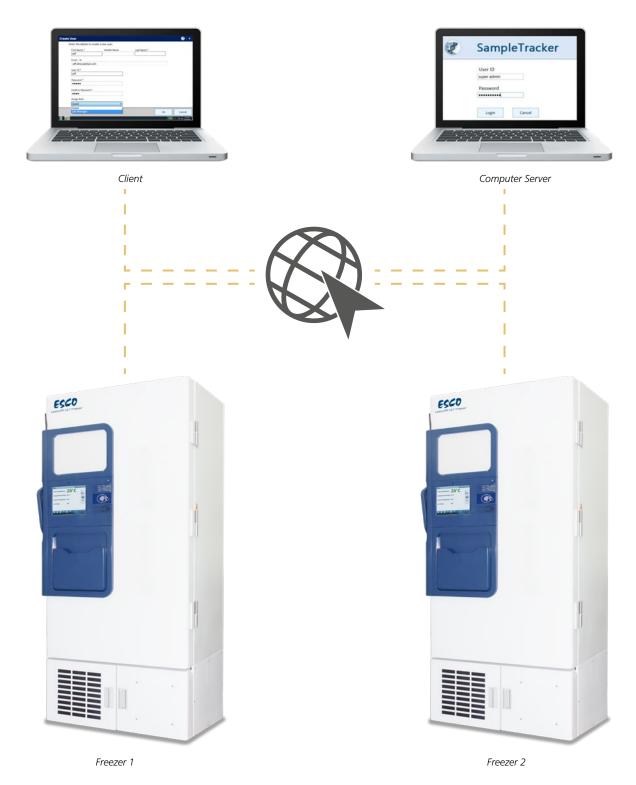


KEY FEATURES

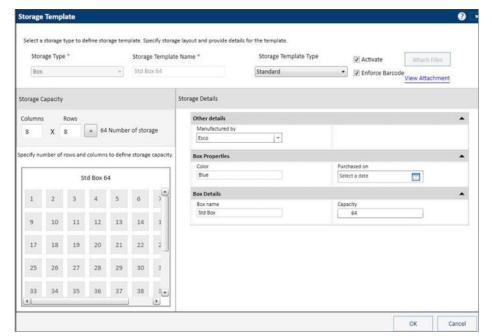


Laboratory Storage Management

- Detail storage definition and configuration
- User able to manage their sample such as save, update and save as another storage
- User will allow to design their own storage template such as capacity and details storage information

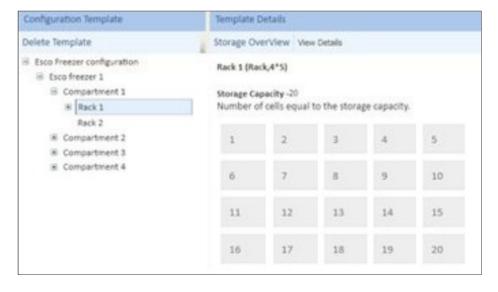

Lab Management

- Detail hierarchies as a configuration template
- navigate through all the configuration templates listed

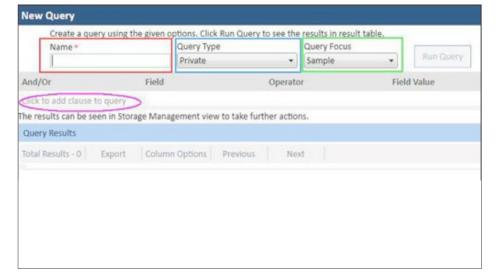

Query Builder

- User will allow to create their own query to search the specific data from database
- User can save the most popular queries and can use these queries in future instead of creating new query every time



User allow to mark their sample that will help them for identifying the individual label in the million samples



SampleTracker / Alto Platinum Controller functions



Login to SampleTracker

At the Alto Platinum controller home screen go to application and select Inventory System

Type in username and password at SampleTracker login screen.

Performing SampleTracker Functions

- CHECK IN: It allows user to Check In the samples again into the system after processing. Only Checked Out samples can be Checked In.
- CHECK OUT: It allows user to take out the samples for processing. Only Checked In samples can be Checked out.
- DISPOSE: Dispose option deletes the selected samples. This option can be used to remove unwanted or degraded samples so that these positions are available for new samples. Any one of Checked In, Checked Out or Moved Out State allows disposal of the samples.



Minimum System Requirements

SampleTracker Server

- Operating system: Windows XP, 7 Professional, 8 and Windows 2003-2012 Server. (English)
- Memory: 8 GB
- Hard disk space: 4 GB
- Prerequisite Software: .NET Framework 4.0 (included) SQL Server 2008 R2 Express (included) or higher

SampleTracker Client

- Operating system: Windows XP, 7 Professional, 8 and Windows 2003 and 2008 Server (English)
- Memory: 8 GB
- Hard disk space: 20 GB
- Prerequisite Software: .NET Framework 4.0 (included)

SampleTracker

After Sales Services



Parts Availability

Whenever service is needed and parts are required, minimizing downtime is a critical objective. Statistical usage analysis helps Esco to predict parts life, permitting Esco to manage logistics and stage proper inventories around the world. The combination of predictive maintenance, historical data and geospecific proximity assures our customers that parts and labor are available whenever service is scheduled through the local sales organization.

Registration, Documentation and Instruction

Quality control at Esco extends from research and development through engineering, manufacturing, shipment, delivery and customer feedback. Esco maintains an aggressive program to encourage warranty card registration by mail, email or online submission so that we know where Esco products are located and how they are being used. Rest assured that all information disclosed from warranty registrations will be kept confidential. All Esco products include unique serial numbers for identification. Documentation for all performance tests is archived and maintained for customer reference.

Online Technical Information

Site preparation instructions are useful before product arrival and installation. Installation and start-up manuals, operation manuals and quick reference guides are available anytime from the Esco resources online.

NSF International Accreditations

The National Sanitation Foundation (NSF) International is an independent, non-profit organization that provides standards development, product certification, auditing, education and risk management for public health and the environment.

In line with Esco's commitment in providing world class services worldwide, Esco has a large contingent of NSF accredited certifiers which makes Esco not only an Excellent Standards COmpany but also an Excellent Service COmpany, which exemplifies Esco's collective quest of being an Eternally Successful COmpany.

The NSF mark is your assurance that the product complies with all the standard requirements, tested by one of the most respected independent certification organizations in existence today. NSF conducts periodic unannounced inspections and product testing to verify that the product continues to comply with the standard. It is valued by consumers, manufacturers, retailers and regulatory agencies worldwide.

References and Links

For more information, you can visit Esco at www.escoglobal.com





ESCO GLOBAL NETWORK

43 LOCATIONS IN 23 COUNTRIES ALL OVER THE WORLD





Biological Safety Cabinets

CO₂ Incubators

Compounding Pharmacy Equipment
Containment / Pharma Products

Ductless Fume Hoods

Lab Animal Research Products

Laboratory Centrifuges

Laboratory Fume Hoods

Laboratory Ovens and Incubators

Laboratory Shakers

Laminar Flow Clean Benches

PCR Cabinets

PCR Thermal Cyclers

Powder Weighing Balance Enclosures

Ultra-low Temperature Freezers

The Esco Group of Companies is a global life sciences tools provider with sales in over 100 countries. The group is active in lab equipment, pharma equipment and medical devices. Manufacturing facilities are located in Asia and Europe. R&D is conducted worldwide spanning the US, Europe and Asia. Sales, service and marketing subsidiaries are located in 12 major markets including the US, UK, Singapore, Japan, China and India. Regional distribution centers are located in the US, UK, and Singapore.

Life Science • Chemical Research • Assisted Reproductive Technology (ART) • Pharmaceutical Equipment • General Equipment





WORLD CLASS, WORLDWIDE,

Esco Micro Pte. Ltd. • 21 Changi South Street 1 • Singapore 486 777 Tel +65 6542 0833 • Fax +65 6542 6920 • mail@escoglobal.com www.escoglobal.com

Esco Technologies, Inc. • 903 Sheehy Drive, Suite F, Horsham, PA 19044, USA Tel +1 215-441-9661 • Fax 484-698-7757 eti.admin@escoglobal.com • www.escolifesciences.us

Esco Global Offices: Bangladesh | Cameroon | China | Denmark | Germany | Hong Kong | India | Indonesia | Italy | Japan | Lithuania | Malaysia | Myanmar | Philippines | Russia | Singapore | South Africa | South Korea | Taiwan | Thailand | UAE | UK | USA | Vietnam







